Ranger[®] 7000 Scales

User Guide

EN

ES

FR

DE

IT

PT

SV

DA

PL

NL

CZ

HU

1. SAFETY INFORMATION

Definition of Signal Warnings and Symbols

WARNING	For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
CAUTION	For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
Attention	For important information about the product.
Note	For useful information about the product.

Warning Symbols

General hazard

Electrical shock

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

Intended Use

Use the instrument exclusively for <weighing/moisture determination/etc.> as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended.



This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

2. INSTALLATION

2.1 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes. Allow sufficient space around the instrument.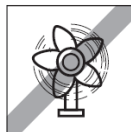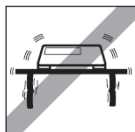

2.2 Connecting Power

AC Power Cord Installation


Connect the AC mains power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

Attention: For optimal weighing performance, allow the scale to warm up for 60 minutes prior to use.

2.3 Connecting the Interface

Use the built-in RS-232 port to connect either to a computer or a printer with a standard (straight-through) serial cable. Or connect using the scale's USB port.

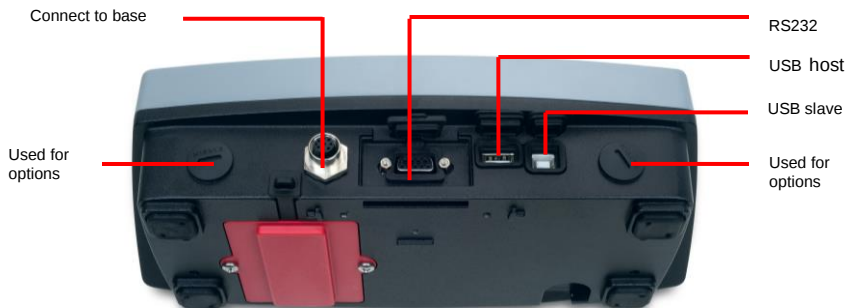Interface connections on back of Terminal.

Thread terminal cable along cable coils on bottom of scale.
Or pass cable through groove near release button.

2.4 Levelling the Instrument

To level the instrument, adjust the feet/leveling wheel so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.

2.5 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

3. OPERATION

3.1 Overview of Display, Home Screen

CONTROLS

Button	Action
	Enter/Exit the library menu
	Switch between available application modes
	Send the measurement data to available communications ports according to current settings.
	Display information about Application Mode, Library, User and Menu
	Enter/Exit the User menu
	Switch the main weighing unit between the available units
	Short Press: Input '2'-'9' To Enter 'A' press 2 times. For lower case 'Z', press 5 times.
	Short Press: Input '0' Long Press: Go to User Login screen
	Short Press: Input '1' Long Press: Switch platform between scale 1 and scale 2
	Short Press: Clear character/string when editing string If no input is active, clear the current active library When there is no value added, pressing this button will switch the value sign between positive and negative.
	Short Press: Input '.', space, '_' To Enter '_' press 3 times.
	Perform Zero operation
	Perform Tare operation When entering the value first and then pressing this button the number input will be set to preset Tare value.

Main Application Screen

MENU & SCREEN NAVIGATION

Press the **Menu**  button to open the menu list.

Press the button below  and  to move down and up the list respectively.

To select the highlighted menu item, press . Press  to move back to previous screen.

3.2 Using the Instrument

Weighing Application

Use this application to determine the weight of items in the selected unit of measure.

Weighing

1. Press the  button until **Weighing** is displayed in the upper left portion of the home screen (this application is the default).
2. Press **Tare** or **Zero** if necessary to begin.
3. Place objects on the pan to display the weight. When stable, the * appears.
4. The resulting value is displayed in the main Weighing Line in the active unit of measure.

4. MAINTENANCE.

4.1 Cleaning

WARNING: Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

4.2 Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website www.ohaus.com to locate the Ohaus office nearest you.

5. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only
- Altitude: Up to 2000 m
- Operating Temperature range: 5 °C to 40 °C.
- Specified Temperature range: 10 °C to 30 °C (R71MHD3/6/15/35 models) -10 °C to 40 °C (R71MD3/6/15/35 models)
- Humidity: maximum relative humidity 80 % for temperatures up to 30 °C decreasing linearly to 50 % relative humidity at 40 °C
- Mains supply voltage fluctuations: up to ± 10 % of the nominal voltage
- Installation category II
- Pollution degree: 2

Specifications:

MODEL	R71MHD3	R71MHD6	R71MHD15	R71MHD35
Capacity	3000 g	6000 g	15000 g	35000 g
Readability d	0.01 g	0.02 g	0.1 g	0.1 g
Approved Readability e	0.1 g	0.2 g	1 g	1 g
Repeatability (std. dev.)	0.01 g	0.02 g	0.1 g	0.1 g
Linearity	± 0.02 g	± 0.04 g	± 0.2 g	± 0.2 g
Power supply	Power Input: 100-240 V~ 0.5 A 50/60 Hz			

MODEL	R71MD3	R71MD6	R71MD15	R71MD35	R71MD60
Capacity	3000 g	6000 g	15000 g	35000 g	60000 g
Readability d	0.05 g	0.1 g	0.2 g	0.5 g	1 g
Approved Readability e	0.5 g	1 g	2 g	5 g	10 g
Repeatability (std. dev.)	0.05 g	0.1 g	0.2 g	0.5 g	1 g
Linearity	± 0.1 g	± 0.2 g	± 0.4 g	± 1 g	± 2 g
Power supply	Power Input: 100-240 V~ 0.5 A 50/60 Hz				

6. COMPLIANCE


Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at www.ohaus.com/ce .
	This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee .

Verified weighing instruments

When the instrument is used in trade or a legally controlled application it must be set up, verified and sealed in accordance with local weights and measures regulations. It is the responsibility of the purchaser to ensure that all pertinent legal requirements are met.

Weighing Instruments verified at the place of manufacture bear the following supplementary metrology marking on the descriptive plate.

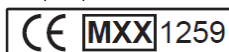Weighing Instruments to be verified in two stages have no supplementary metrology marking on the descriptive plate. The second stage of conformity assessment must be carried out by the applicable weights and measures authorities. If national regulations limit the validity period of the verification, the user of the weighing instrument must strictly observe the re-verification period and inform the weights and measures authorities

As verification requirements vary by jurisdiction, the purchaser should contact their local weights and measures office if they are not familiar with the requirements

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P/N 30283470 A © 2015 Ohaus Corporation, all rights reserved.
Ohaus Corporation